Guardian 5000 / 7000 Quick Start Guide

EN

ES

FR

DE

IT

PT

SE

NL

DA

FI

NO

PL

CZ

HU

BASIC FUNCTIONS

Quick Press

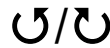
Long Press

Press & Hold

Rotate

<p>I/O – Standby Switch</p>	<p>Turn unit on (Press switch)</p>	
	<p>Turn unit off (Press switch)</p>	
<p>Temp Knob Functions</p>	<p>Change temp settings (Rotate temp knob)</p>	
	<p>Turn heater on (Long press temp knob until unit beeps & green light appears)</p>	
	<p>Change temperature settings, while heating (Rotate temp knob, the setting will blink, then quick press knob)</p>	<p>(If setting is idle for 6 sec, setting will return to prior value)</p>
	<p>Turn heater off (Long press temp knob until unit beeps & green light fades)</p>	
	<p>Enter settings menu <i>[heater & stirrer must be off]</i> (Press & hold temp knob until “MENU” appears)</p>	
<p>Speed Knob Functions</p>	<p>Change speed settings (Rotate speed knob)</p>	
	<p>Turn stirrer on (Long press speed knob until unit beeps & green light appears)</p>	
	<p>Change speed settings, while stirring (Rotate speed knob, the setting will blink, then quick press knob)</p>	<p>(If setting is idle for 6 sec, setting will return to prior value)</p>
	<p>Turn stirrer off (Long press speed knob until unit beeps & green light fades)</p>	
	<p>Enter timer settings - Guardian 7000 only <i>[heater & stirrer must be off]</i> (Press & hold speed knob until “HH:MM” appears)</p>	

MENU (Not applicable for Stirrers [e-G51ST07C])

To access the settings menu, the heater and stirrer must be off. Then press and hold the Temp/Menu knob until “MENU” appears on the screen. (The heater will not turn on unless the knob is released before “MENU” appears.)

To navigate the settings menu, rotate the knob then quick press the knob to choose the selection. Once in the selection, rotate the knob to change the setting and quick press the knob to confirm.

To return to the top level of the settings menu, rotate the knob to “BACK” then quick press the knob. To return to the main operating screen from the top level of the settings menu, rotate the knob to “EXIT” then quick press the knob.

To exit the menu at any time, flip the standby switch off and back on. The unit will return to the main operating screen and will be ready for normal use.

Menu Features:

1. “TLIM” - SmartHeat™ (Guardian 7000 units only)

SmartHeat™ allows the user to change the maximum temperature limit of the heater.

2. “PROG” – Programs (Guardian 7000 units only)

The Programs menu allows the user to load, store, delete, and run a series of heater and stirrer functions from a programmable text file on a USB storage device.

3. “RAMP” – SmartRate™ (Guardian 7000 units only)

SmartRate™ allows the user to change the rate at which the heater and/or stirrer increases temperature or speed respectively.

4. “SAFE” – SmartPresence™ & SmartLink™ (Guardian 7000 units only)

SmartPresence™ is an optional safety feature designed to automatically turn the heater off if the unit is unattended for a user-selected time limit.

SmartLink™ is an optional safety feature designed to automatically turn the heater off if the user travels out of range of the unit for a user-selected time limit.

Note: SmartLink™ requires the optional Ohaus Wireless Dongle accessory.

5. “CAL” – Single Point Calibration

Single Point Calibration (SPC) improves the accuracy of the heater at user-selected temperature points. SPC can be used during plate-controlled and probe-controlled heating.

6. “SYS” – System Settings

- i. Enabling / Disabling the Beeper
- ii. Changing the Timer Start Setting (Guardian 7000 units only)
- iii. Changing the Power Recovery Setting
- iv. Resetting to Factory Default Settings

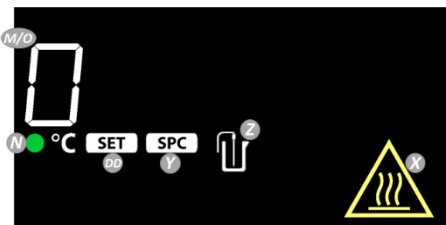
DISPLAY (Guardian 7000)

DISPLAY (Guardian 5000)

DISPLAY (Guardian 5000 Hotplate 7 x 7)

DISPLAY (Guardian 5000 Stirrer 7 x 7)

- L. Heater Indicator Bar:** Illuminates when the heater is $\geq 40^{\circ}\text{C}$.
- M. Heater Temperature:** Switches to external probe temperature when the probe is plugged in and Z is illuminated.
- N. Heater Indicator:** Illuminates when heater is running
- O. Heat Setting**
- P. Stir Speed**
- Q. Stirrer Indicator:** Illuminates when stirrer is running
- R. Speed Setting**
- S. Temperature Priority Icon**
- T. Timer:** Hours : Minutes / Minutes : Seconds
- U. Program Icon:** Illuminates when a program is running.
- V. SmartPresence™ Icon**
- W. Bluetooth® Icon:** Illuminates when *SmartLink™* is enabled.
- X. Hot Top Caution Indicator:** Illuminates when the heater is $\geq 40^{\circ}\text{C}$.
- Y. Single Point Calibration Icon**
- Z. External Probe Icon**
- AA. SmartHeat™ Icon**
- BB. SmartRate™ Heater Icons:**
- Turtle –slower ramp rate (more precise)
 - Rabbit –faster ramp rate
- CC. SmartRate™ Stirrer Icons:**
- Turtle –slower ramp rate (more precise)
 - Rabbit –faster ramp rate
- DD. Heat Setting Indicator:** Switches Heater Temperature to Heat Setting when illuminated.
- EE. Speed Setting Indicator:** Illuminates until stirrer reaches the Speed Setting.

SAFETY INFORMATION

Safety notes are marked with signal words and warning symbols. These show safety issues and warnings. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results.

WARNING	For a hazardous situation with medium risk, possibly resulting in severe injuries or death if not avoided.
CAUTION	For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or minor or medium injuries if not avoided.
ATTENTION	For important information about the product. May lead to equipment damage if not avoided.
NOTE	For useful information about the product.

Warning Symbols

General
Hazard

Caution, hot
surface

Electrical
shock hazard

SAFETY PRECAUTIONS

WARNING: DO NOT use the Hotplate-Stirrer in explosive atmospheres or with materials that could cause a hazardous environment from processing. Keep in mind the material flash point relative to the target temperature that has been set. Also, the user should be aware that the protection provided by the equipment may be impaired if used with accessories not provided by the manufacturer.

Always operate unit on a level surface for best performance and maximum safety.

DO NOT lift unit by the top plate.

CAUTION! To avoid electrical shock, completely cut off power to the unit by disconnecting the power cord from the wall outlet. Disconnect unit from the power supply prior to maintenance and servicing.

Spills should be removed promptly after the unit has cooled down.

DO NOT immerse the unit for cleaning. Alkalis spills, hydrofluoric acid or phosphoric acid spills may damage the unit and lead to thermal failure.

CAUTION! The top plate can reach 500°C.

DO NOT touch the heated surface. Use caution at all times. Keep the unit away from explosive vapors and clear of papers, drapery, and other flammable materials. Keep the power cord away from the heater plate.

CAUTION! The rear panel of the 10x10-120V unit runs hot to the touch. Avoid contact during operation. Allow unit to cool before touching the rear panel.

DO NOT operate the unit at high temperatures without a vessel/sample on the top plate.

DO NOT operate the unit if it shows sign of electrical or mechanical damage.

WARNING! Units are NOT explosion proof. Use caution when heating volatile materials.

Earth Ground – Protective Conductor Terminal. Protective earthing of the equipment is achieved via connection of the provided power cord to a compatible grounded power outlet.

Alternating Current

COMPLIANCE

Compliance to the following standards is indicated by the corresponding mark on the product.

Mark	Standard
	OHAUS Corporation declares that the Guardian series hotplates, stirrers, and hotplate-stirrers comply with directives 2011/65/EU, (EU) 2015/863, 2014/30/EU, 2014/35/EU, and standards EN 50581, EN 61010-2-010, EN 61010-2-051, EN 61326-1. The full text of the EU declaration of conformity is available at the following internet address: www.ohaus.com/ce .
	This product complies with directive 2012/19/EU. Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. For disposal instructions in Europe, refer to www.ohaus.com/weee .
	EN 61326-1
	CAN/CSA C22 261010-1, CAN/CSA C22 261010-2-010, CAN/CSA C22 261010-2-051 UL 61010-1, UL 61010-2-010, UL 61010-2-051

Global Notice

Warning: This is a Class A product. In a domestic environment this product may cause radio interference in which case the user may be required to take adequate measures.

Canada Notice

This Class A digital apparatus complies with Canadian ICES-003.

FCC Notice

NOTE: This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

Changes or modifications not expressly approved by Ohaus Corporation could void the user's authority to operate the equipment.

TROUBLESHOOTING

The following table lists common problems and possible causes and remedies. If the problem persists, contact OHAUS or your authorized dealer.

Error*	Cause of Error	How to Fix
Unit fails to power on	Missing or blown fuse	Add or replace fuse as necessary.
E1	Plate RTD open	Not fixable by user, please contact Ohaus.
E2	Plate RTD short	Not fixable by user, please contact Ohaus.
E3	No stirring motion / cannot reach speed	Not fixable by user, please contact Ohaus.
E4	Probe RTD open (Removing the probe while the unit is heating)	Switch unit to standby, then return to normal operating mode.
E5	Probe RTD short (Malfunctioning probe)	Switch unit to standby, remove the probe from the unit, then return to normal operating mode.
E6	A/D lock error	Not fixable by user, please contact Ohaus.
E7	User Probe Error (Plugging the probe into the unit while it is heating)	Switch unit to standby, then return to normal operating mode.
E8	Plate over temperature	Not fixable by user, please contact Ohaus.
E9	Plate under temperature	Not fixable by user, please contact Ohaus.
E10	Triac fault	Not fixable by user, please contact Ohaus.

***Note:** Error code instances will stop equipment operation by default.

INSTRUCTION MANUAL

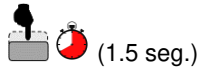
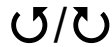
For the full instruction manual, please scan the QR code or click the links below: You can also visit <http://www.ohaus.com/support> to download user manual.

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SAP P/N 30500607A P/N 786501-00 © 2020 Ohaus Corporation, all rights reserved