

SE572

Installation, Maintenance & Operation Instructions

DESCRIPTION

Deck mounted emergency eyewash. Supplied with 15mm stay-open full flow ball valve; activated by stainless steel push handle; twin spray outlets with flow restrictor, and flow off dust covers. Unit swivels 90° for easy access to sink when not required, and mounts on counter tops up to 1-1/4" thick. **Manufactured in Australia to comply with AS 4775 - 2007.**

SPECIFICATIONS

SPRAY HEADS: (2) Aerated Stainless Steel with Dust Cover.

VALVE: 15mm Stainless Steel Ball Valve

ACTIVATOR: Stainless Steel Push Handle.

INLET: 1/2" BSP

OPTIONS INSTALLATION

At desired location, drill a 25mm (1") diameter hole through countertop (1-1/4" thick max.) Be sure the hole is drilled as straight as possible. Attach eyewash to countertop using supplied hardware, and connect a water supply to unit that is a 12mm minimum uninterrupted source of potable water (use pipe sealant on connection). Failure to use potable water can result in emergency units producing impure or contaminated water causing possible further injury.



TESTING PROGRAM

The AS 4775 - 2007 states that all safety emergency equipment shall be activated on a weekly basis to flush the line and verify proper operation. All units furnishes a testing tag (SE Tag) . On this tag the date of the inspection and the inspectors initials should be noted.

NOTE: This unit should be connected to an uninterrupted source of potable water, with a minimum flowing pressure of 30 Kpa and a maximum static pressure of 125 PSI.

OPERATION

Be sure unit is above sink or drain before activating. The eyewash is activated by pushing the push handle back 90°. Once the unit is activated, water will discharge from the spray outlets until the push handle is manually pulled back to the off position.

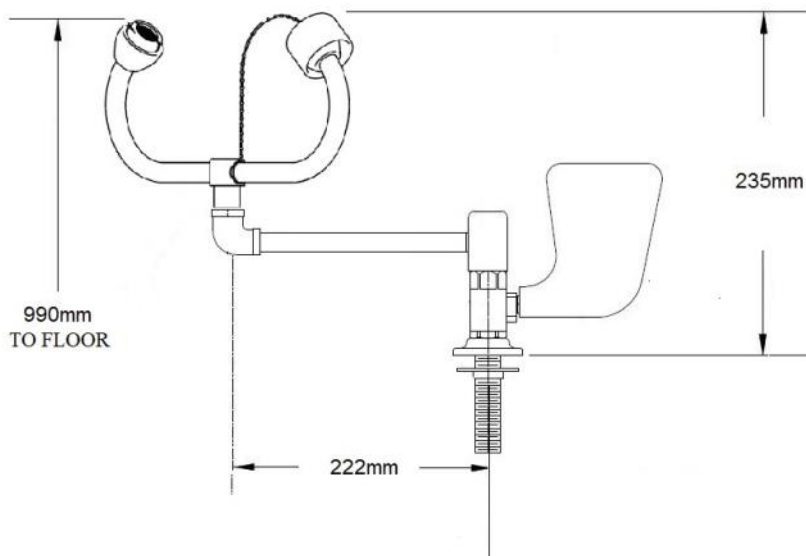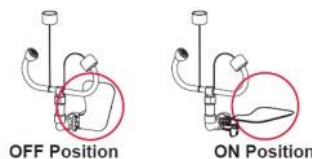

EYE WASH FLOW RATE (LPM)

WARNING : Use only genuine Speakman parts when replacing or repairing components. When ordering quote correct part numbers listed.

SPEAKMAN®
SAFETY PRODUCTS

SE 572

Swivel Bench Mount Eye Wash

Installation, Maintenance & Operation Instructions

MAINTENANCE

The valve and the eyewash aerators are the only portions of the unit that may require maintenance. Should you need parts to repair this unit, please reference the parts listing below for correct part numbers.

MAINTENANCE SCHEDULE


WEEKLY	Activate eyewash unit using push handle Allow water to flow for 30 seconds to flush line. Ensure water flows evenly from both outlets. Ensure water is clean.
YEARLY	Ensure unit complies with AS 4775-2007

- 3 Part No. RPG05 1177
Aerator Screen & Cap Assembly




Sold as a packet of 6.

- 1 Part No: SA 01003

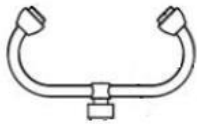

Complete Bull Horn Assembly

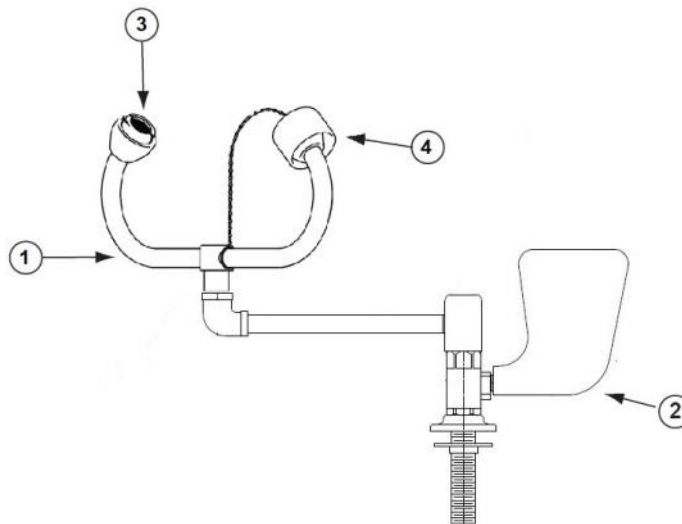

SPARE PARTS

Item	Description	Part No
1.	Complete Aerated Bullhorn Assy.	SA 01003
2.	15mm Ball Valve & Push Handle Ass	SE 910V
3.	Aerator Screen & Cap Assembly	RPG05 1177
4.	Red Dust Cap & Chain	SA 01008

WARNING : Use only genuine Speakman parts when replacing or repairing components. When ordering quote correct part numbers listed.

